- 
- (EN) Bird Rinse**
- 
- 
- (FR) Flacon pour rinçage de volaille**
- 
- 
- (DE) Geflügelspülwasser**
- 
- 
- (IT) Bird Rinse**
- 
- 
- (ES) Enjuague para Aves**
- 
- 
- (NL) Bird Rinse**
- 
- 
- (SV) Fågel-rinse**
- 
- 
- (DA) Bird Rinse**
- 
- 
- (NO) Fugleprøve**
- 
- 
- (FI) Lintunäytteiden huuhteluaine**
- 
- 
- (PT) Bird Rinse**
- 
- 
- (EL) Έκπλυμα Πτηνών**
- 
- 
- (PL) Płukanka do ptaków**
- 
- 
- (HU) Bird Rinse szárnyasmosó**
- 
- 
- (CS) Láhev Bird Rinse**
- 
- 
- (RO) Bird Rinse**
- 
- 
- (RU) Раствор для промывания птицы**
- 
- 
- (TR) Bird Rinse**
- 
- 
- (JA) 食鳥洗淨液**
- 
- 
- (ZH) Bird Rinse**
- 
- 
- (TH) Bird Rinse**
- 
- 
- (KO) 조류용 린스**

# Product Instructions

## Bird Rinse

### PRODUCT DESCRIPTION AND INTENDED USE

3M™ Bird Rinse bottles come pre-filled with 400 mL of either Butterfield's Buffer or Buffered Peptone Water Broth. 3M Bird Rinse will save time when used with the 3M™ Carcass kits. The wide bottle opening makes broth dispensing easy and the seal provides assurance that the product has not been previously opened.

**Table 1: Product Description and Storage Conditions**

Catalog Number	Product Description	Volume	Storage Symbol
BRBPW400	3M™ Bird Rinse with Buffered Peptone Water Broth	400 mL	20-25°C
BRBFD400	3M™ Bird Rinse with Butterfield's Buffer	400 mL	20-25°C

3M Food Safety is certified to ISO (International Organization for Standardization) 9001 for design and manufacturing

### SAFETY

The user should read, understand, and follow all safety information in the instructions for the 3M Bird Rinses. Retain the safety instructions for future reference.

**⚠ CAUTION** Indicates a hazardous situation, which, if not avoided, could result in minor or moderate injury and/or property damage.

<b>⚠ CAUTION</b>
------------------

- 3M has not documented 3M Bird Rinse diluents for use in industries other than food. For example, 3M has not documented 3M Bird Rinse for testing water, pharmaceuticals, or cosmetics.
- 3M Bird Rinse diluents have not been tested with all possible food products, food processes, testing protocols or with all possible strains of bacteria.
- Do not use 3M Bird Rinse in the diagnosis of conditions in humans or animals.
- The user must train its personnel in proper testing techniques, for example: Good Laboratory Practices<sup>(3)</sup>, ISO 7218<sup>(4)</sup> or ISO 17025<sup>(5)</sup>.
- For information on documentation of product performance, visit our website at [www.3M.com/foodsafety](http://www.3M.com/foodsafety) or contact your local 3M representative or distributor.
- Consult the Safety Data Sheet for additional information.

### USER RESPONSIBILITY

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at [www.3M.com/foodsafety](http://www.3M.com/foodsafety), or contact your local 3M representative or distributor for more information.

When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results. The food sample itself may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples with the appropriate matrices and microbial challenges to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any 3M Food Safety product do not constitute a guarantee of the quality of the matrices or processes tested.

### LIMITATION OF WARRANTIES / LIMITED REMEDY

EXCEPT AS EXPRESSLY STATED IN A LIMITED WARRANTY SECTION OF INDIVIDUAL PRODUCT PACKAGING, 3M DISCLAIMS ALL EXPRESS AND IMPLIED WARRANTIES, INCLUDING BUT NOT LIMITED TO, ANY WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE. If any 3M Food Safety Product is defective, 3M or its authorized distributor will, at its option, replace or refund the purchase price of the product. These are your exclusive remedies. You must promptly notify 3M within sixty days of discovery of any suspected defects in a product and return it to 3M. Please call Customer Service (1-800-328-1671 in the U.S.) or your official 3M Food Safety representative for a Returned Goods Authorization.

### LIMITATION OF 3M LIABILITY

3M WILL NOT BE LIABLE FOR ANY LOSS OR DAMAGES, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING BUT NOT LIMITED TO LOST PROFITS. In no event shall 3M's liability under any legal theory exceed the purchase price of the product alleged to be defective.

### STORAGE

Upon receipt, store unopened 3M Bird Rinse at the indicated temperatures (See Table 1).

## **⚠ DISPOSAL**

After use, 3M Bird Rinse may contain microorganisms that may be a potential biohazard. Follow current industry standards or local regulations for disposal.

## **INSTRUCTIONS FOR USE**

Follow all instructions carefully. Failure to do so may lead to inaccurate results.

Wear appropriate protective apparel and follow standard good laboratory practices GLP<sup>1</sup>.

1. Unscrew and remove cap from 3M Bird Rinse bottle.
2. Put on gloves.
3. To remove seal, grasp the tab and peel back from bottle edge, continuing to pull the seal away from the bottle until removed.
4. Tear open a 3M™ Bird Rinse Bag (BRB3500) or equivalent.
5. Pour the 3M Bird Rinse diluent into the bag containing the carcass cavity.
6. Follow your defined protocol for bird carcass rinsing<sup>(1, 2)</sup>.
7. After rinsing, pour the sample rinse fluid back into the bottle, close cap.

## **REFERENCES**

1. ISO 17604 AMD 1:2009. Microbiology of food and animal feeding stuffs - Carcass sampling for microbiological analysis.
2. USDA FSIS MLG 3.01. Quantitative Analysis of Bacteria in Foods as Sanitary Indicators, Section 3.5.3 Whole Bird Rinse.
3. U.S. Food and Drug Administration. Code of Federal Regulations, Title 21, Part 58. Good Laboratory Practice for Nonclinical Laboratory Practice Studies.
4. ISO 7218. Microbiology of food and animal feeding stuffs – General requirements and guidance for microbiological examinations.
5. ISO/IEC 17025. General requirements for the competence of testing and calibration laboratories.

Refer to the current versions of the standard methods listed above.

## **EXPLANATION OF SYMBOLS**

Consult product instructions.

The lot in a box and the hourglass symbols are symbols that represent lot number and expiration date. The hourglass is followed by a year, month, and day which represents the expiration date (year, month, and day: YYYY-MM-DD). The entire line after the hourglass represents the lot number (YYYY-MM AZ).

Store between given temperatures.

## 3M Food Safety

### 3M United States

3M CenterBldg. 275-5W-05  
St. Paul, MN 55144-1000  
USA  
1-800-328-6553

### 3M Canada

Post Office Box 5757  
London, Ontario N6A 4T1  
Canada  
1-800-563-2921

### 3M Europe and MEA

3M Deutschland GmbH  
Carl-Schurz - Strasse 1  
D41453 Neuss/Germany  
+49-2131-14-3000

### 3M Latin America

3M Center  
Bldg. 275-5W-05  
St. Paul, MN 55144-1000  
USA  
1-954-340-8263

### 3M Asia Pacific

No 1, Yishun Avenue 7  
Singapore, 768923  
65-64508869

### 3M Japan

3M Health Care Limited  
6-7-29, Kita-Shinagawa  
Shinagawa-ku, Tokyo  
141-8684 Japan  
81-570-011-321

### 3M Australia

Bldg A, 1 Rivett Road  
North Ryde, NSW 2113  
Australia  
61 1300 363 878

### 3M Health Care

2510 Conway Ave  
St. Paul, MN 55144 USA  
[www.3M.com/foodsafety](http://www.3M.com/foodsafety)

© 2014, 3M. All rights reserved.  
3M and Petrifilm are trademarks of 3M. Used under license in Canada.  
34-8715-5735-0

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1. ISO 17604 AMD 1:2009. Microbiology of food and animal feeding stuffs - Carcass sampling for microbiological analysis.
2. USDA FSIS MLG 3.01. Quantitative Analysis of Bacteria in Foods as Sanitary Indicators, Section 3.5.3 Whole Bird Rinse.
3. U.S. Food and Drug Administration. Code of Federal Regulations, Title 21, Part 58. Good Laboratory Practice for Nonclinical Laboratory Studies.
4. ISO 7218. Microbiology of food and animal feeding stuffs – General requirements and guidance for microbiological examinations.
5. ISO/IEC 17025. General requirements for the competence of testing and calibration laboratories.

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## REFERINȚE

1. ISO 17604 AMD 1:2009. Microbiology of food and animal feeding stuffs - Carcass sampling for microbiological analysis.
2. USDA FSIS MLG 3.01. Quantitative Analysis of Bacteria in Foods as Sanitary Indicators, Section 3.5.3 Whole Bird Rinse.
3. U.S. Food and Drug Administration. Code of Federal Regulations, Title 21, Part 58. Good Laboratory Practice for Nonclinical Laboratory Practice Studies.
4. ISO 7218. Microbiology of food and animal feeding stuffs - General requirements and guidance for microbiological examinations.
5. ISO/IEC 17025. General requirements for the competence of testing and calibration laboratories.

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© 2015, 3M. All rights reserved.