Clean-Trace™

Quick Start Guide

- Ⓔ **EN** Hygiene Monitoring and Management System Quick Start Guide
- Ⓕ **FR** Guide de démarrage rapide du contrôle d'hygiène et système de management
- Ⓖ **DE** Hygiene Monitoring und Management System – Kurzanleitung
- Ⓙ **IT** Guida rapida per il Sistema di gestione e di verifica dell'igiene
- Ⓚ **ES** Guía de inicio rápido del Sistema de Monitoreo y Gestión de Higiene
- Ⓝ **NL** Snelstartgids Hygiëne Monitoring en Management Systeem
- Ⓢ **SV** Snabbstartguide för Hygiene Monitoring and Management System
- Ⓓ **DA** Kvikstartguide til Hygiejne Monitorerings System
- Ⓝ **NO** Oppstartsveiledning for system for hygienekontroll
- Ⓕ **FI** Hygienian valvonta- ja hallintajärjestelmän pikaopas
- Ⓕ **PT** Guia de Início Rápido do Monitoramento de Higiene e Sistema de Gestão
- Ⓔ **EL** Σύστημα Ελέγχου και Διαχείρισης της Υγιεινής - Σύντομος Οδηγός Εκκίνησης
- Ⓕ **PL** Skrócona instrukcja obsługi systemu monitorowania higieny i zarządzania higieną
- Ⓕ **HU** Higiéniamonitorozó és -menedzsment rendszer - Rövid beüzemelési útmutató
- Ⓕ **CS** Systém monitorování a řízení hygieny – stručný návod
- Ⓕ **RO** Ghid rapid de utilizare al sistemului de monitorizare și administrare a igienei
- Ⓕ **RU** Краткое руководство по началу работы с системой мониторинга чистоты
- Ⓕ **TR** Hijyen Mönitörizasyon ve Yönetim Sistemi Hızlı Başlangıç Kılavuzu
- Ⓕ **JA** 衛生モニタリング/マネジメントシステム クイックスタートガイド
- Ⓕ **ZH** 卫生监控系统快速入门指南 (Chinese, Simplified)
- Ⓕ **ZH** 衛生監控管理系統快速入門指引 (Chinese, Traditional)
- Ⓕ **TH** คู่มือการใช้งานฉบับย่อสำหรับระบบการจัดการการตรวจสอบประสิทธิภาพการทำงานสะอาด
- Ⓕ **KO** 환경위생 검사 및 관리 시스템 빠른 시작 안내서
- Ⓕ **ID** Panduan Mulai Cepat Hygiene Monitoring and Management System

3M™ Clean-Trace™ Hygiene Monitoring and Management System Quick Start Guide (EN)

Consisting of 3M™ Clean-Trace™ Luminometer and 3M™ Clean-Trace™ Hygiene Management Software

PACKAGE CONTENTS

1. 3M™ Clean-Trace™ Luminometer
2. 3M™ Clean-Trace™ Hygiene Management Software (Stored on 3M Clean-Trace Luminometer)
3. 3M™ Clean-Trace™ Luminometer Base
4. Power supply with interchangeable country plugs
5. USB 2.0 mini-A communications cable
6. 3M Clean-Trace Hygiene Monitoring and Management System Quick Start Guide
7. 3M™ Clean-Trace™ Hygiene Monitoring and Management System Important Information booklet
8. Certificate of Conformity

1. Read and follow all Safety Information in the 3M Clean-Trace Hygiene Monitoring and Management System Important Information booklet
2. 3M™ Clean-Trace™ Hygiene Monitoring and Management System User Manual

 Available at www.3M.com/CleanTraceLM1manual

3. Charge the 3M Clean-Trace Luminometer for at least 2 hours before initial use.

Select appropriate power supply adapter; connect cord to power supply, connect power supply to 3M Clean-Trace Luminometer


4. Confirm your computer meets the minimum requirements

- Microsoft® Windows® 7 (32-bit or 64-bit), or Microsoft® Windows® 8 (32-bit or 64-bit) or Microsoft® Windows® 8.1 (32- or 64-bit)
- 2.0 GHz Intel Pentium 4 or similar processor
- 1 GB RAM
- 10 GB of free disk space
- USB 2.0 compatible

IMPORTANT! Complete steps 1-4 before installing and launching 3M Clean-Trace Hygiene Management Software

5. Connect USB cable to computer and power on 3M Clean-Trace Luminometer by depressing the Power button (⏻) for 3 seconds.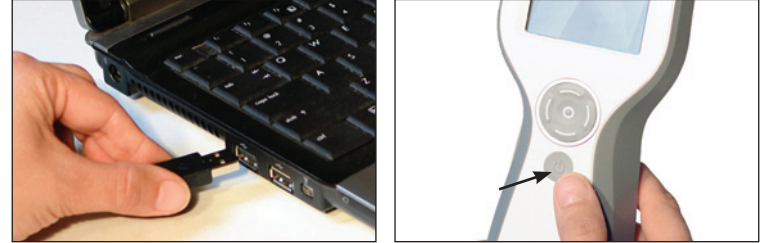

6. Install the 3M Clean-Trace Hygiene Management Software by following the installation instructions located at www.3M.com/CleanTraceLM1manual in the 3M Clean-Trace Hygiene Monitoring and Management System User Manual

Disclaimer: 3M disclaims all express and implied warranties including warranties of merchantability or fitness for a particular use. If product is defective, the exclusive remedy is, at 3M's option, replacement, or refund.

Except where prohibited by law, 3M will not be liable for any further loss or damage arising from use of this product.

User Responsibility: Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.3M.com/foodsafety, or contact your local 3M representative or distributor for more information.

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3M and Clean-Trace are trademarks of
All other trademarks are the property of
©3M 2016. All rights reserved.
3M Health Care Limited