-  (EN) Transfer Pipette
-  (FR) Pipette Jetable
-  (DE) Transfer Pipette
-  (IT) Pipetta di trasferimento
-  (ES) Pipeta de Transferencia
-  (NL) Transferpipet
-  (SV) Transfer Pipette
-  (DA) Pipette
-  (NO) Overføringspipette
-  (PT) Pipeta de Transferência
-  (EL) Πιπέτα Δειγματοληψίας
-  (PL) Pipeta
-  (RU) Трансферная пипетка
-  (TR) Pipetör
-  (JA) トランスファーピペット
-  (ZH) 移液管
-  (TH) ปิเปตต์ถ่ายตัวอย่าง
-  (KO) 1회용 스포이드
-  (ID) Pipet Transfer
-  (AR) ماصة نقل

Product Instructions

Transfer Pipette

Product Description and Intended Use

The 3M™ Transfer Pipette is intended for use with the 3M™ Allergen Protein Rapid Kits to transfer clean-in-place (CIP) final rinse water to the 3M™ Extraction Buffer, or any extracted sample to the 3M™ Allergen Protein Lateral Flow Device (LFD). The 3M Transfer Pipette package contains 100 transfer pipettes.

Safety

The user should read, understand and follow all safety information in the product instructions for the 3M Transfer Pipette. Retain the safety instructions for future reference.

⚠ WARNING: Indicates a hazardous situation, which, if not avoided, could result in death or serious injury and/or property damage.

⚠ NOTICE: Indicates a potentially hazardous situation, which, if not avoided, could result in property damage.

⚠ WARNING

To reduce the risks of inaccurate results:

- 3M has not documented the use of this product in industries other than food or beverage. For example, 3M has not documented this product for testing pharmaceutical, cosmetic, clinical or veterinary samples.
- 3M Transfer Pipette should be used to load 0.1 mL of the sample diluted in the 3M™ Extraction Buffer onto the 3M Allergen Protein LFD.
- 3M Transfer Pipette may be used to transfer twice 0.1 mL of CIP final rinse water but its re-use with other samples should be verified.
- Store the 3M Transfer Pipette as indicated on the package and in the product instructions.
- Always use the 3M Transfer Pipette by the expiration date.
- Always use the 3M Transfer Pipette at 20-30°C.

3M Transfer Pipette is intended for use in the Food and Beverage industries by trained personnel. For information on documentation of product performance, visit our website at www.3M.com/foodsafety or contact your local 3M representative or distributor.

User Responsibility

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.3M.com/foodsafety, or contact your local 3M representative or distributor for more information.

As with all test methods used for food analysis the test matrix can influence the results. When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results. The food sample itself may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any 3M Food Safety product do not constitute a guarantee of the quality of the matrices or processes tested.

Limitation of Warranties/Limited Remedy

EXCEPT AS EXPRESSLY STATED IN A LIMITED WARRANTY SECTION OF INDIVIDUAL PRODUCT PACKAGING, 3M DISCLAIMS ALL EXPRESS AND IMPLIED WARRANTIES, INCLUDING BUT NOT LIMITED TO, ANY WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE. If any 3M Food Safety Product is defective, 3M or its authorized distributor will, at its option, replace or refund the purchase price of the product. These are your exclusive remedies. You must promptly notify 3M within sixty days of discovery of any suspected defects in a product and return it to 3M. Please call Customer Service (1-800-328-1671 in the U.S.) or your official 3M Food Safety representative for a Returned Goods Authorization.

Limitation of 3M Liability

3M WILL NOT BE LIABLE FOR ANY LOSS OR DAMAGES, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING BUT NOT LIMITED TO LOST PROFITS. In no event shall 3M's liability under any legal theory exceed the purchase price of the product alleged to be defective.

Storage and Disposal

Store 3M Transfer Pipette at 2 to 30°C.

3M Transfer Pipette should not be used past the expiration date. Expiration date and lot number are noted on the outside label of the box.

Dispose according to current local/regional/industry standards and regulations.

Instructions for Use

Follow all instructions carefully. Failure to do so may lead to inaccurate results. The 3M Transfer Pipette should be used at ambient temperature (20-30°C).

The description of the 3M Transfer Pipette is shown in Figure A.

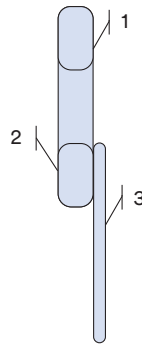
Figure A. Parts of the 3M Transfer Pipette. (1) Upper bulb, (2) Overflow reservoir, (3) Sample tube.

After combining the sample with the 3M Extraction Buffer per the product instructions in the 3M Allergen Protein Rapid Kits, use the 3M Transfer Pipette to transfer 0.1 mL of the prepared sample to the 3M Allergen Protein LFD, as shown in Figure B.

For CIP rinse water only, if using the 3M Transfer Pipette to transfer 0.2 mL of CIP final rinse water sample to the 3M Extraction Buffer, using the same transfer pipette, repeat the steps to transfer 0.1 mL once more for a final transferred volume of 0.2 mL. Use a new 3M Transfer Pipette to transfer 0.1 mL of the prepared sample to the 3M Allergen Protein LFD as shown in Figure B.

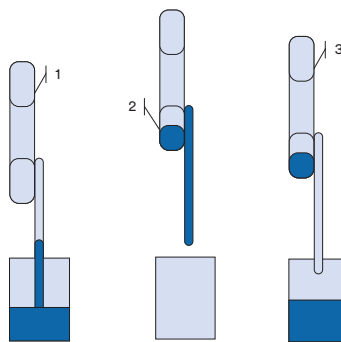
Figure B. Procedure to collect 0.1 mL utilizing the 3M Transfer Pipette. (1) Squeeze upper bulb to completely fill sample tube. (2) Excess sample will flow into overflow reservoir. (3) Squeeze upper bulb to dispense liquid from sample tube. Overflow liquid will remain in overflow reservoir.

Explanation of Symbols

www.3M.com/foodsafety/symbols

3M Food Safety

3M United States

3M Center
Bldg. 275-5W-05
St. Paul, MN 55144-1000
USA
1-800-328-6553

3M Canada

Post Office Box 5757
London, Ontario N6A 4T1
Canada
1-800-563-2921

3M Latin America

3M Center
Bldg. 275-5W-05
St. Paul, MN 55144-1000
USA
1-954-340-8263

3M Europe and MEA

3M Deutschland GmbH
Carl-Schurz-Strasse 1
D41453 Neuss/Germany
+49-2131-14-3000

3M United Kingdom PLC

Morley Street, Loughborough
Leicestershire
LE11 1EP
United Kingdom
+(44) 1509 611 611

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3M Health Care

2510 Conway Ave
St. Paul, MN 55144 USA
www.3M.com/foodsafety

© 2018, 3M. All rights reserved.
3M and Petrifilm are trademarks of 3M. Used under license in Canada.
34-8723-4135-8

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3M Österreich GmbH
3M Health Care Limited