

# 3M™ Harvest RC Chromatographic Clarifier


Single-stage chromatographic  
purification for recombinant  
protein therapeutic  
manufacturing



Clarification using 3M™ Harvest RC Chromatographic Clarifier provides many benefits including fewer processing steps, higher yield and lower production time and costs.

This single-stage single-use chromatographic clarification solution is the next generation in harvest and clarification technology. It is designed as an efficient option for harvesting and clarification of modern cell cultures.


**3M™ Harvest RC Chromatographic Clarifier improves process economics with:**

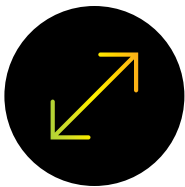


Scalability from bench to  
commercial production




Predictable  
performance

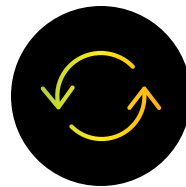

High  
product recovery

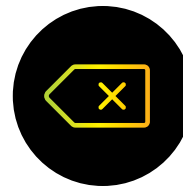

Whole cell  
removal and minimize  
cell shear

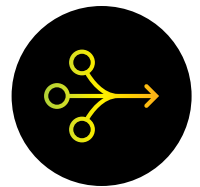

Streamlined  
process and  
smaller footprint

## Innovative design and performance

3M™ Harvest RC Chromatographic Clarifier encapsulates innovative synthetic fibrous anion exchange (AEX) chromatography media and a 0.2 µm polyether sulfone (PES) membrane. This enables a single-stage clarification process of low- to high-density cell culture (>40 million cells per mL) with high recovery, and high fidelity of soluble and insoluble contaminant separation.

Cells are bound inside the media by electrostatic charge interaction with the AEX chromatographic fibres. This results in the efficient retention of large and small particulates without developing a surface cake layer. The media can also remove soluble impurities which results in cleaner effluent than centrifugation or depth filtration.



*Before and after using 3M™ Harvest RC Chromatographic Clarifier: turbidity reduction in a single stage*

### Key features and benefits:

- ▶ Simplification of high cell density cell culture fluid clarification unit operations
- ▶ Optimized for PCV high density CHO cell culture (5-8% PCV)
- ▶ Replacement of primary, secondary, and guard membrane clarification stages
- ▶ Typical product recovery of 95+% (capsules)
- ▶ Synthetic chromatographic harvest media with chemically defined extractables
- ▶ Predictable scaling from discovery to manufacturing
- ▶ Lower total cost of manufacturing compared to centrifugation and depth filtration
- ▶ Lower consumption of buffer and water compared to depth filtration
- ▶ Capsules fit into laboratory to manufacturing scale workflows




# 3M™ Harvest RC Chromatographic Clarifier – Single step chromatographic clarification encapsulated solution


The innovative synthetic fibrous AEX chromatographic clarification media enables a single-stage clarification process of low- to high-density CHO cell culture (> 40 million cells per mL) with high product recovery, and high fidelity of soluble and insoluble contaminant separation.

Downstream of the fibrous chromatographic clarification media is the 0.2 µm PES membrane which distributes the flow across the AEX media bed and enables protection of downstream sterilizing grade membrane filter. Also, 0.2 µm PES membrane enables simple process endpoint measurement using pressure reading.


## Expanding fibrous media platform

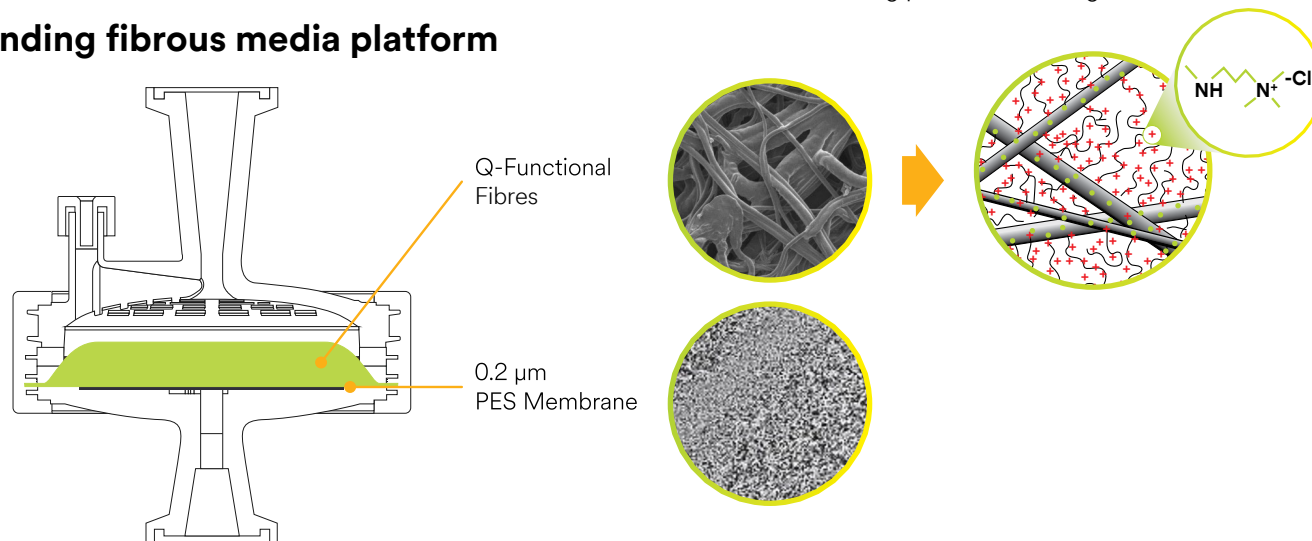

## Cell harvest simplified:

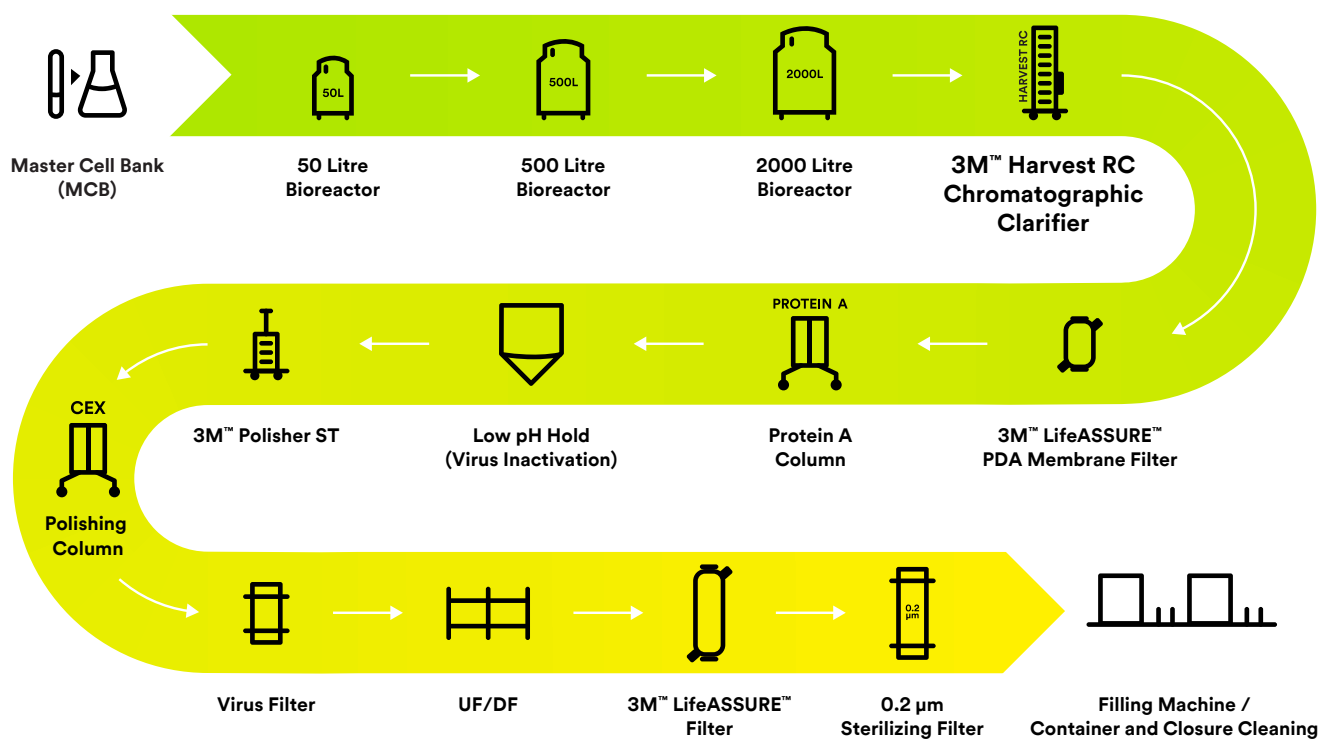

# The next generation in harvest and clarification technology

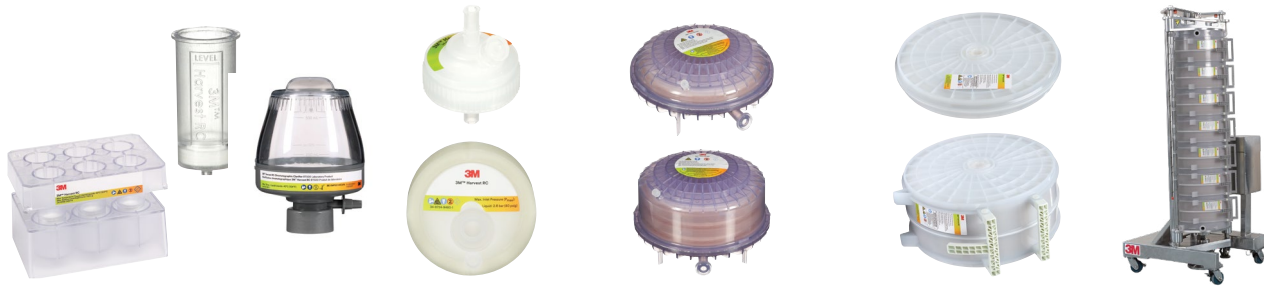

Research & Development

Scale-up

Manufacturing

## 3M™ Harvest RC Chromatographic Clarifier product family: laboratory, pilot, and production capsules



### Efficient workflows. Excellent clarity.

With 3M™ Harvest RC Chromatographic Clarifier, BT500, it's just 10 minutes to a high-quality sample for downstream processing.\* No centrifuge is needed, and the easy-to-use, bottle-top vacuum clarifier enables high product recoveries and DNA reduction (<500 ppb). Together, these properties let you streamline your upstream and downstream processes before they start, which eliminates complexity and improves workflow.

\*Based on 5-8% PCV culture




## Performance data

All performance data below is based on capsules. While it is representative of typical performance, results may vary depending on the format(s) used.


### mAb product recovery

3M™ Harvest RC Chromatographic Clarifier is a single stage chromatography solution that effectively clarifies Chinese Hamster Ovary (CHO) harvest cell culture fluid (HCCF) across a wide range of cell densities, packed cell volumes (PCV), and turbidities.

3M™ Harvest RC Chromatographic Clarifier capsules consistently provide >95% mAb product recovery for high cell density cultures from the laboratory to the manufacturing scale.

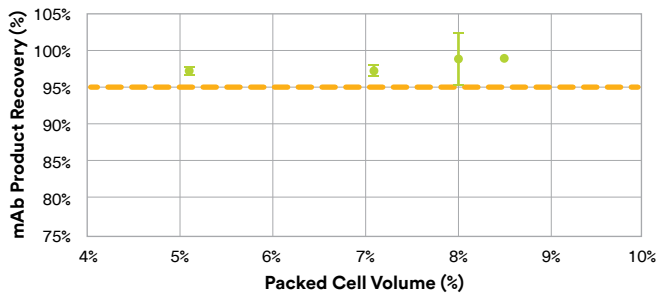


**Figure 1A:** mAb product recovery in clarification process at different packed cell volumes (N = 1 – 4)

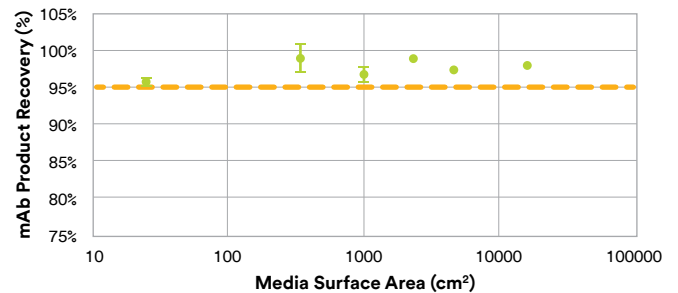

**Figure 1B:** mAb product recovery in clarification process at different media surface areas (N = 1 – 3)

### Turbidity reduction

3M™ Harvest RC Chromatographic Clarifier provides consistent separation of cells, cell debris, and DNA from the target protein. Clarified cell culture fluid (CCCF) has low turbidity, typically <15 NTU. Additionally, consistently low acidified turbidity of CCCF indicates significant reduction of DNA in the clarified material. Low acidified CCCF turbidity is a measure of the amount of DNA present in the cell culture fluid. (Koehler et al. Biotechnology Progress. 2019;35:e2882)

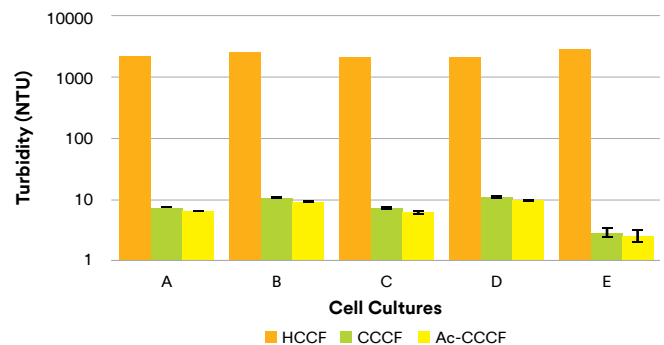


**Figure 2:** Turbidity Reduction by 3M™ Harvest RC capsules (N = 3 – 6). A – E are different CHO cell cultures at 5 – 8 % PCV.

### Scalability

3M™ Harvest RC Chromatographic Clarifier capsules scale linearly across laboratory, pilot, and manufacturing scales.

Fibrous chromatographic clarification assures scalable performance from discovery to manufacturing scales. Performance is consistent from laboratory capsules (BC4 and BC25), scale-up capsules (BC340 and BC1020), to production capsules (BC2300 and BC16000) within  $\pm 20\%$  of BC25 throughput.

Throughputs of 3M™ Harvest RC Chromatographic Clarifier capsules are scaled by area based on packed cell volume.

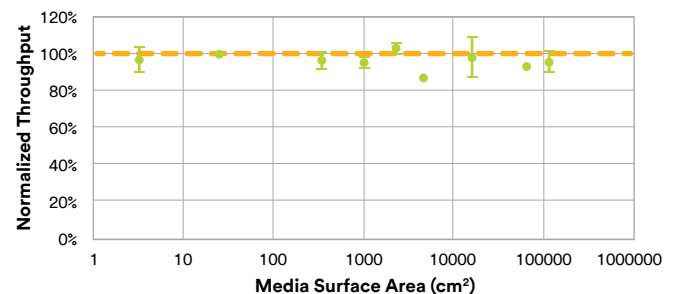


**Figure 3:** Scalability from laboratory to scale-up and production capsules (N = 1 – 5, 6 cell cultures)

## Performance data (continued)

### Cell loading capacity


3M™ Harvest RC Chromatographic Clarifier utilizes advanced Q functionalized fibrous chromatography media to achieve single-stage clarification, enabling predictable and consistent cell loading capacity for CHO cell culture fluid for a wide range of packed cell volumes.

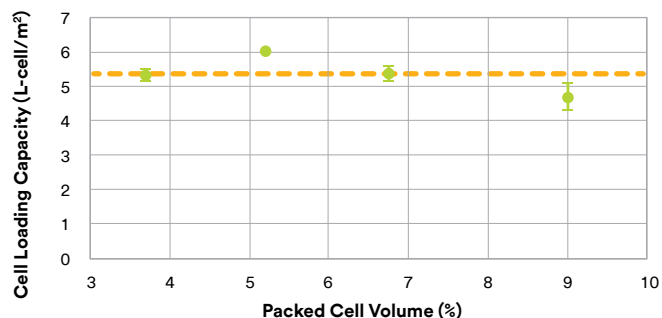

**Figure 4:** Cell loading capacity of 3M™ Harvest RC capsules for CHO harvested cell culture fluid at different packed cell volumes (N = 2 - 3)

### Cell shear


The low-pressure chromatographic clarification relies on charge rather than size or density. This results in minimal cell shear compared to conventional depth filtration processes even at medium and high cell densities. Cell shear was evaluated by lactate dehydrogenase (LDH) assay (Sigma-Aldrich 11644793001).

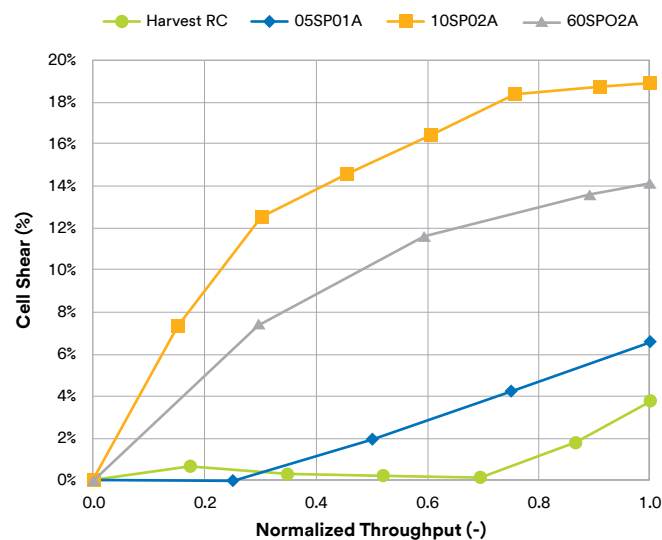

**Figure 5:** Minimal cell shear of 3M™ Harvest RC during clarification of 8% PCV CHO cell culture at 100 LMH.

### Robust sterile filter protection

Due to the highly effective chromatographic reduction of soluble and insoluble contaminants, 3M™ Harvest RC Chromatographic Clarifier enables efficient clarification, and is capable of effective protection of final sterilizing grade membrane filter down to 0.1 μm pore size.

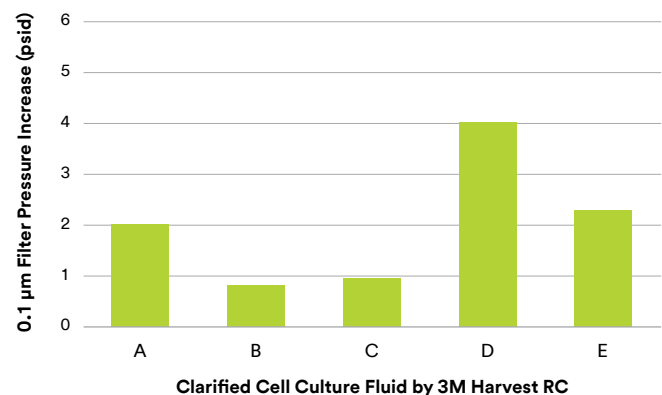

**Figure 6:** 0.1 μm sterile filter pressure increase at 500 L/m<sup>2</sup>. A – E are clarified fluids of CHO harvested cell culture fluids at 8%PCV by 3M™ Harvest RC capsules.

# Product specifications



Product Name	BC4	BC25 Luer	BC25 Sanitary	BC340	BC1020	BC2300	BC16000
<b>Model Name</b>	EMP201HRC2FA	EMP301HRC2FA	EMP303HRC2FA	EMP513HRC2FA	EMP533HRC2FA	EMP710HRC2FA	EMP770HRC2FA
<b>Global Part Number</b>	70-0203-5331-7	70-0203-5332-5	70-0203-5333-3	70-0203-5335-8	70-0203-5336-6	70-0203-5337-4	70-0203-5339-0
<b>EMA Part Number</b>	7100241969	7100236863	7100236867	7100236865	7100236866	7100236845	7100236846
<b>Height x Diameter</b>	5.9 cm x 4.3 cm (2.3 in x 1.7 in)	5.3 cm x 7.7 cm (2.1 in x 3.0 in)	8.6 cm x 7.7 cm (3.4 in x 3.0 in)	10.4 cm x 24.1 cm (4.1 in x 9.5 in)	15.2 cm x 24.1 cm (6.0 in x 9.5 in)	5.7 cm x 45.2 cm (2.2 in x 17.8 in)	20.3 cm x 45.2 cm (8.0 in x 17.8 in)
<b>Dry Weight</b>	14.3 g	69.2 g	75.8 g	1.1 kg	1.6 kg	3.4 kg	9.8 kg
<b>Media Surface Area</b>	3.2 cm <sup>2</sup>	25 cm <sup>2</sup>	25 cm <sup>2</sup>	340 cm <sup>2</sup>	1020 cm <sup>2</sup>	2300 cm <sup>2</sup>	1.61 m <sup>2</sup>
<b>Cell Culture Volume Range (5-8% PCV)<sup>1</sup></b>	20 - 32 mL	150 - 250 mL	150 - 250 mL	2 - 3.4 L	6 - 10 L	14 - 23 L	100 - 160 L
<b>Weight Wet Post Blow Down</b>	17.2 g	81.2 g	88.1 g	1.2 kg	2.1 kg	4.4 kg	16.3 kg
<b>Fill Volume<sup>2</sup></b>	5.6 mL	27.6 mL	28.2 mL	0.66 L	1.7 L	3.3 L	16.3 L
<b>Hold up Volume Post Blow Down<sup>3</sup></b>	3.0 mL	12.0 mL	12.3 mL	0.16 L	0.47 L	1.1 L	6.5 L
<b>Capsule Material</b>	Polypropylene	Polypropylene, Glass Filled Polypropylene	Polypropylene, Glass Filled Polypropylene	Polysulfone, Polypropylene, Glass Filled Polypropylene, Thermoplastic Elastomer, Fluorocarbon	Polysulfone, Polypropylene, Glass Filled Polypropylene, Thermoplastic Elastomer, Fluorocarbon	Polycarbonate, Polypropylene, Glass Filled Polypropylene, Thermoplastic Elastomer, Silicone	Polycarbonate, Polypropylene, Glass Filled Polypropylene, Thermoplastic Elastomer, Silicone
<b>Inlet / Outlet Connections</b>	Luer-Lok	Luer-Lok	Sanitary	Sanitary	Sanitary	Sanitary	Sanitary
<b>Maximum Inlet Pressure<sup>4</sup></b>	3.4 bar	2.8 bar	2.8 bar	3.1 bar	3.1 bar	3.4 bar	3.4 bar
<b>Maximum Differential Pressure</b>	2.4 bar	2.4 bar	2.4 bar	2.4 bar	2.4 bar	2.4 bar	2.4 bar
<b>Maximum Temperature</b>	40 °C (104 °F)	40 °C (104 °F)	40 °C (104 °F)	40 °C (104 °F)	40 °C (104 °F)	40 °C (104 °F)	40 °C (104 °F)
<b>Required Preconditioning Flush Volume<sup>5</sup></b>	8 mL	62.5 mL	62.5 mL	0.85 L	2.55 L	5.8 L	40.3 L
<b>Recommended Use Flow Rate</b>	0.53 mL/min	4.2 mL/min	4.2 mL/min	57 mL/min	170 mL/min	0.38 L/min	2.68 L/min
<b>Storage Conditions</b>	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging
<b>Shelf Life</b>	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature
<b>cGMP Compliant</b>	No	No	No	Yes	Yes	Yes	Yes
<b>GLP Compliant</b>	Yes	Yes	Yes	No	No	No	No



Product Name	WP6	CT15	BT500
<b>Model Name</b>	EMP006HRC2FA	EMP015HRC2FA	EMP051HRC2FA
<b>Global Part Number</b>	70-0203-5328-3	70-0203-5329-1	70-0203-6542-8
<b>EMA Part Number</b>	7100242700	7100245350	7100302871
<b>Height x Diameter</b>	12.8 cm x 8.5 cm x 8.8 cm (5.0 in x 3.4 in x 3.5 in)	2.9 cm x 6.1 cm (1.2 in x 2.4 in)	16.1 cm x 11.1 cm (6.3 in x 4.4 in)
<b>Dry Weight</b>	Plate (w/ Media): 100 g Collector Plate: 110 g	10 g	250 g
<b>Cell Culture Volume Range (5-8% PCV)<sup>1</sup></b>	15 mL per well	15 mL	500 mL
<b>Fill Volume<sup>2</sup></b>	15 mL per well	15 mL	500 mL
<b>Capsule Material</b>	Polycarbonate	Polycarbonate	Polycarbonate
<b>Maximum Temperature</b>	40 °C (104 °F)	40 °C (104 °F)	40 °C (104 °F)
<b>Maximum Relative Centrifugal Force</b>	750 x g	750 x g	N/A
<b>Recommended Relative Centrifugal Force</b>	400 x g	400 x g	N/A
<b>Recommended Spin Time</b>	10 minutes	10 minutes	10 minutes
<b>Maximum Vacuum Pressure</b>	N/A	N/A	15 psi
<b>Recommended Vacuum Pressure</b>	N/A	N/A	7.5 psi
<b>Storage Conditions</b>	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging	Controlled indoor temperatures: 0-30 °C (32-86 °F) in original sealed packaging
<b>Shelf Life</b>	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature	Up to 2 years from the date of manufacture @30 °C maximum storage temperature
<b>cGMP Compliant</b>	No	No	No
<b>GLP Compliant</b>	Yes	Yes	Yes

## Footnotes:

- Cell Culture Volume Range is the estimation for CHO cell culture fluid at 5 - 8% packed cell volume
- Fill Volume is defined as the volume of liquid that is required to fill the capsule.
- Post Blow-Down Hold-Up Volume is defined as the volume of the residual liquid after air/gas blow down
- Do not use this product for continuous service with compressed gasses. The use of compressed gas is permissible for post-use integrity testing and blow down purposes.
- A Preconditioning Flush is required for the product to be compliant with USP Biological Reactivity Tests, including USP <87> and <88> Class VI. Refer to Installation and Operation Instructions for complete instructions on how to perform the preconditioning flush.

For more information about the **3M™ Harvest RC Chromatographic Clarifier**, contact your local sales or application engineering representative by calling **1-800-243-6894**, option 4, or visiting us at **3M.com/Bioprocessing**

**Intended Use:** 3M™ Harvest RC products are intended for use in biopharmaceutical processing applications of aqueous based pharmaceuticals (drugs) and vaccines in accordance with the product instructions and specifications, and cGMP requirements (for BC340, BC1020, BC2300 and BC16000) or GLP requirements (for CT15, WP6, BC4 and BC25), where applicable.

Since there are many factors that can affect a product's use, the customer and user remain responsible for determining whether the 3M product is suitable and appropriate for the user's specific application, including user conducting an appropriate risk assessment and evaluating the 3M product in user's application.

**Restrictions on Use:**

*For CT15, WP6, BC4 and BC25:* For laboratory use only. Not intended for use with materials that will be used on humans or animals.

For all sizes: 3M advises against the use of these 3M products in any application other than the stated intended use(s), since other applications have not been evaluated by 3M and may result in an unsafe or unintended condition. Do not use in any manner whereby the 3M product, or any leachable from the 3M product, may become part of or remains in a medical device that is regulated by any agency, and/or globally exemplary agencies, including but not limited to: a) FDA, b) European Medical Device Regulation (MDR), c) Japan Pharmaceuticals and Medical Devices Agency (PMDA) or in applications involving permanent implantation into the body; Life-sustaining medical applications; Applications requiring food contact compliance.

**Technical Information:** The technical information, guidance, and other statements contained in this document or otherwise provided by 3M are based upon records, tests, or experience that 3M believes to be reliable, but the accuracy, completeness, and representative nature of such information is not guaranteed. Such information is intended for people with knowledge and technical skills sufficient to assess and apply their own informed judgment to the information. No license under any 3M or third party intellectual property rights is granted or implied with this information.

**Product Selection and Use:** Many factors beyond 3M's control and uniquely within user's knowledge and control can affect the use and performance of a 3M product in a particular application. As a result, end-user is solely responsible for evaluating the product and determining whether it is appropriate and suitable for end-user's application, including completing a risk assessment that considers the product leachable characteristics and its impact on drug safety, conducting a workplace hazard assessment and reviewing all applicable regulations and standards. Failure to properly evaluate, select, and use a 3M product and appropriate safety products, or to meet all applicable safety regulations, may result in injury, sickness, death, and/or harm to property.

**Warranty, Limited Remedy, and Disclaimer:** Unless a different warranty is expressly identified on the applicable 3M product literature or packaging (in which case such express warranty governs), 3M warrants that each 3M product meets the applicable 3M product specification at the time 3M ships the product. 3M MAKES NO OTHER WARRANTIES OR CONDITIONS, EXPRESS OR IMPLIED, INCLUDING, BUT NOT LIMITED TO, ANY IMPLIED WARRANTY OR CONDITION OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR ARISING OUT OF A COURSE OF DEALING, CUSTOM, OR USAGE OF TRADE. If a 3M product does not conform to this warranty, then the sole and exclusive remedy is, at 3M's option, replacement of the 3M product or refund of the purchase price.

**Limitation of Liability:** Except for the limited remedy stated above, and except to the extent prohibited by law, 3M will not be liable for any loss or damage arising from or related to the 3M product, whether direct, indirect, special, incidental, or consequential (including, but not limited to, lost profits or business opportunity), regardless of the legal or equitable theory asserted, including, but not limited to, warranty, contract, negligence, or strict liability.



**3M Separation and Purification Sciences Division**

300 Tartan Drive  
London, ON N5V 4M9  
Canada

1-800-364-3577  
3M.ca

3M, 3M Science. Applied to life. and LifeASSURE are trademarks of 3M. Used under license in Canada. © 2023, 3M. All rights reserved. All other trademarks are the property of their respective owners. 2310-27051 E