

Cleaning LifeASSURE™ BA Series Membrane Filters with Caustic Solutions

SAFETY INFORMATION

Read, understand, and follow all safety information contained in these instructions and the instructions provided with the original filtration system, prior to installation and use. Retain for future reference.

EXPLANATION OF SIGNAL WORD CONSEQUENCES	
WARNING:	Indicates a potentially hazardous situation, which, if not avoided, could result in death or serious injury and/or property damage.
CAUTION:	Indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury and/or property damage.

WARNING:
To reduce the risk associated with system burst related injuries: <ul style="list-style-type: none"> • Do not use if fluid pressure exceeds rating described on the pressure vessel dataplate. • Do not use with fluids at temperatures exceeding the rating described on the pressure vessel dataplate. • Do not use for continuous service with compressed gases.

CAUTION:
To reduce the risk associated with exposure to contaminants: <ul style="list-style-type: none"> • Always use appropriate personal protective equipment (PPE) when installing or servicing the filtration system. • Ensure that all system pressure has been relieved prior to opening the system to atmosphere. • To reduce the risk associated with eye, skin, and respiratory and digestive tract injuries from chemical cleaners/sanitizers during system maintenance: <ul style="list-style-type: none"> ○ Do not get chemical cleaners/sanitizers in eyes, on skin, or on clothing. Do not ingest or inhale. ○ Wear appropriate PPE including eye and face protection, protective gloves, and an appropriate NIOSH-approved filter mask.

Introduction

In malt or fruit based beverage applications where pre-filtration is used prior to the final membrane filters, the final membrane filters tend to plug by gradual constriction of the pores with colloidal substances, not by the collection of particulate or microorganisms on the membrane surface. The colloidal substances are composed of a variety of materials that include beta-glucans, proteins and tannins. These components are inherent in the malt or fruit, or are generated in upstream processes such as fermentation and the enzymatic breakdown of carbohydrates and other compounds.

Regular flushing of the filter with warm water can reduce this build up (refer to 3M Purification Inc. Technical Brief 70-0202-5340-0). However, flushing with a dilute caustic solution can even further reduce colloidal build-up. This procedure has been found to work best with beer or malt based beverages. As an example, in the filtration of beer, laboratory studies show a 25% minimum filter service life increase when using a warm water flush in conjunction with a mild alkaline flush as opposed to the warm water flush alone. For best results, it is essential to clean the filters before the process flow differential pressure exceeds 10 psid (~700 mbar).



The recommended alkaline solution is 2-3% NaOH at a maximum of 140°F (60°C). Always use filtered water to make the solution.

Procedure:

Warning: Exercise caution when working with caustic solutions. Appropriate apparel, gloves, and face/eye protection (safety glasses, goggles, or full mask) must be worn at all times.

1. At the end of a standard filtration cycle, flush the filter with warm 131 °F (55 °C) filtered water as described in 3M Purification Technical Brief 70-0202-5340-0. Flush until filtrate appears clear.
2. Drain the housing and close the inlet and outlet housing valves.
3. Open the upstream housing bleed valve.
4. Slowly open the caustic solution valve at the inlet to the housing.
5. Slowly allow the caustic solution to fill the housing.
6. As the caustic solution emerges, close the upstream housing bleed valve and slowly open the downstream housing valve.
7. Flush the caustic solution through the system for 15 minutes at a flow rate at least as fast as the beverage itself (typically 1-2 gpm [4-8 lpm] per 10" filter cartridge). Faster flows will access more pores but do not exceed 3 gpm (12 lpm) per 10" filter cartridge.
8. Flush housing to neutralize any residual alkaline solution in the housing with filtered ambient temperature 0.5% to 1% citric acid solution.
9. Integrity test the filter(s) before next use.

Important Notice:

The test results described in this literature are accurate to the best of our knowledge. A variety of factors, however, can affect the performance of the product(s) in a particular application, some of which are uniquely within your knowledge and control. **INFORMATION IS SUPPLIED UPON THE CONDITION THAT THE PERSONS RECEIVING THE SAME WILL MAKE THEIR OWN DETERMINATION AS TO ITS SUITABILITY FOR THEIR USE. IN NO EVENT WILL 3M PURIFICATION INC. BE RESPONSIBLE FOR DAMAGES OF ANY NATURE WHATSOEVER RESULTING FROM THE USE OF OR RELIANCE UPON INFORMATION.**

It is your responsibility to determine if additional testing or information is required and if the 3M Purification Inc. product is fit for a particular purpose and suitable in your specific application.

3M PURIFICATION INC. MAKES NO REPRESENTATIONS OR WARRANTIES, EITHER EXPRESS OR IMPLIED INCLUDING WITHOUT LIMITATION ANY WARRANTIES OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE OR OF ANY OTHER NATURE HEREUNDER WITH RESPECT TO INFORMATION OR THE PRODUCT TO WHICH INFORMATION REFERS.

Limitation of Liability:

3M Purification Inc. will not be liable for any loss or damage arising from the use of the product(s), whether direct, indirect, special, incidental, or consequential, regardless of the legal theory asserted, including warranty, contract, negligence or strict liability. Some states and countries do not allow the exclusion or limitation of incidental or consequential damages, so the above limitation or exclusion may not apply to you.



3M Purification Inc.
400 Research Parkway
Meriden, CT 06450
U.S.A.
(800) 243-6894
(203) 237-5541
Fax (230) 630-4530
www.3Mpurification.com

3M is a trademark of 3M Company.
LifeASSURE is a trademark of 3M Company used under license.
© 2010 3M Company. All rights reserved.
Please recycle. Printed in the U.S.A.
70-0202-6594-1 REV 1110