

3M™ Petrifilm™ *Salmonella* Express Plate

Presumptive Positive *Salmonella* Colors

Presumptive positive *Salmonella* species are red to brown colonies with a yellow zone or associated gas bubble, or both.

70-2009-9523-4



Note: This color chart serves as a sample of colony colors. Not all colony colors are represented below.



3M™ Petrifilm™ *Salmonella* Express Plate

Non-*Salmonella* Colors

Non-*Salmonella* species are blue, blue to green, green and/or black colonies with or without a yellow zone and/or associated gas bubble.



Note: This color chart serves as a sample of colony colors. Not all colony colors are represented below.

