- (EN) ***Salmonella* Enrichment Base**
- (FR) ***Salmonella* Base pour Enrichissement**
- (DE) ***Salmonella* Anreicherung - Basis**
- (IT) **Base di arricchimento per *Salmonella***
- (ES) **Enriquecimiento Base para *Salmonella***
- (NL) ***Salmonella* Verrijkingsmedium**
- (SV) ***Salmonella* anrikningsbuljong**
- (DA) ***Salmonella* Opformerings Base**
- (NO) ***Salmonella* oppformeringsbuljong**
- (FI) ***Salmonella* Rikastusalusta**
- (PT) **Base para Enriquecimento de *Salmonella***
- (EL) **Σαλμονέλλα, Βάση Εμπλουτισμού**
- (PL) ***Salmonella* Baza wzbogacająca**
- (HU) ***Salmonella* dúsító bázis**
- (CS) **Obohacené médium pro stanovení *Salmonelly***
- (RO) **Bază de îmbogățire pentru *Salmonella***
- (RU) **Среда обогащения для бактерий рода *Salmonella***
- (TR) ***Salmonella* Zenginleştirme Bazi**
- (JA) **サルモネラエンリッチメントベース**
- (ZH) **沙门氏菌基础增菌培养基**
- (TH) **อาหารเลี้ยงเชื้อสำหรับเชื้อซาลโมเนลล่า**
- (KO) **살모넬라 증균 베이스**

Product Instructions

Salmonella Enrichment Base

DESCRIPTION

Used for the selective enrichment of *Salmonella* species in conjunction with the 3M™ Petrifilm™ *Salmonella* Express (SALX) System. Supplement required.

SAFETY

The user should read, understand, and follow all safety information in the instructions for the 3M *Salmonella* Enrichment Base. Retain the safety instructions for future reference.

⚠ WARNING Indicates a hazardous situation, which, if not avoided, could result in death or serious injury and/or property damage.

NOTICE Indicates a potentially hazardous situation which, if not avoided, could result in property damage.

⚠ WARNING

To reduce the risks associated with false-negative results leading to release of contaminated food or beverage product:

- Always follow the proper enrichment scheme (time, temperature, and dilution).
- Always use care when pipetting or transferring samples.
- Follow the protocol and perform the tests exactly as stated in the product instructions for use.

To reduce the risks associated with false negative or false positive results leading to the release of contaminated product or the destruction of good product:

- Always follow product instructions for preparation, handling, and storage.

To reduce the risks associated with exposure to chemicals and biohazards:

- Dispose of enriched samples according to all applicable government regulations and applicable laboratory procedures.
- Read, understand, and follow the safety information in the 3M Petrifilm *Salmonella* Express System Instructions for Use and the associated MSDS.
- Always follow standard good laboratory safety practices (GLP)¹, including proper containment procedures, wearing appropriate protective apparel and eye protection while handling testing materials and test samples.

To reduce the risks associated with environmental contamination:

- Follow current industry standards and local regulations for disposal of contaminated waste.

NOTICE: To reduce false-positive results due to cross contamination that result in re-testing or the rejection of food or beverage product:

- Always wear gloves.
- Use a new pipette tip for each sample transfer.

USER RESPONSIBILITY

Users are responsible for familiarizing themselves with product instructions and information. Visit our website at www.3M.com/foodsafety, or contact your local 3M representative or distributor for more information.

When selecting a test method, it is important to recognize that external factors such as sampling methods, testing protocols, sample preparation, handling, and laboratory technique may influence results.

It is the user's responsibility in selecting any test method or product to evaluate a sufficient number of samples with the appropriate matrices and microbial challenges to satisfy the user that the chosen test method meets the user's criteria.

It is also the user's responsibility to determine that any test methods and results meet its customers' and suppliers' requirements.

As with any test method, results obtained from use of any 3M Food Safety product do not constitute a guarantee of the quality of the matrices or processes tested.

LIMITATION OF WARRANTIES / LIMITED REMEDY

EXCEPT AS EXPRESSLY STATED IN A LIMITED WARRANTY SECTION OF INDIVIDUAL PRODUCT PACKAGING, 3M DISCLAIMS ALL EXPRESS AND IMPLIED WARRANTIES, INCLUDING BUT NOT LIMITED TO, ANY WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR USE. If any 3M Food Safety Product is defective, 3M or its authorized distributor will, at its option, replace or refund the purchase price of the product. These are your exclusive remedies. You must promptly notify 3M within sixty days of discovery of any suspected defects in a product and return it to 3M. Please call Customer Service (1-800-328-1671 in the U.S.) or your official 3M Food Safety representative for a Returned Goods Authorization.

LIMITATION OF 3M LIABILITY

3M WILL NOT BE LIABLE FOR ANY LOSS OR DAMAGES, WHETHER DIRECT, INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES, INCLUDING BUT NOT LIMITED TO LOST PROFITS. In no event shall 3M's liability under any legal theory exceed the purchase price of the product alleged to be defective.

STORAGE AND DISPOSAL

Storage: At ambient temperature (15-25°C). Hygroscopic. Keep container tightly closed. Protect from light, heat, and moisture.

Disposal: After use, enriched samples may contain microorganisms that may be a potential biohazard.

Follow current industry standards and local regulations for disposal of contaminated waste. Consult the Material Safety Data Sheet for additional information.

INSTRUCTIONS FOR USE

Follow all instructions carefully. Failure to do so may lead to inaccurate results.

Wear appropriate protective apparel and follow standard good laboratory safety practices (GLP)¹.

Preparation of Medium: Suspend 37.0 g of the powder in 1.0 L of purified water. Mix thoroughly. If necessary, heat to completely dissolve powder. Dispense. Autoclave at 121°C for 15 minutes.

Supplement required: Refer to 3M™ *Salmonella* Enrichment Supplement (CAT: SESUP001) instructions for use for further information.

Final pH 7.0 ± 0.2 at 25°C.

Composition: (typical formula g/L):

Nutrient Mix 36.5

Selective mix 0.5 g

REFERENCES

1. U.S. Food and Drug Administration. Code of Federal Regulations, Title 21, Part 58. Good Laboratory Practice for Nonclinical Laboratory Studies.

EXPLANATION OF SYMBOLS

Consult instructions for use

The lot in a box and the hourglass symbols are symbols that represent lot number and expiration date. The hourglass is followed by a year, month, and day which represents the expiration date (year, month, and day: YYYY- MM-DD). The entire line after the LOT represents the lot number.

Storage conditions

3M Food Safety

3M United States

3M Center
Bldg. 275-5W-05
St. Paul, MN 55144-1000
USA
1-800-328-6553

3M Canada

Post Office Box 5757
London, Ontario N6A 4T1
Canada
1-800-563-2921

3M Europe and MEA

3M Deutschland GmbH
Carl-Schurz - Strasse 1
D41453 Neuss/Germany
+49-2131-14-3000

3M Latin America

3M Center
Bldg. 275-5W-05
St. Paul, MN 55144-1000
USA
1-954-340-8263

3M Asia Pacific

No 1, Yishun Avenue 7
Singapore, 768923
65-64508869

3M Japan

3M Health Care Limited
33-1, Tamagawadai 2-chrome
Setagaya-ku, Tokyo
158-8583, Japan
81-570-011-321

3M Australia

Bldg A, 1 Rivett Road
North Ryde, NSW 2113
Australia
61 1300 363 878

3M Health Care

2510 Conway Ave
St. Paul, MN 55144 USA
www.3M.com/foodsafety

© 2013, 3M. All rights reserved.
3M and Petrifilm are trademarks of 3M. Used under license in Canada.
34-8711-1225-5

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